

System/Product	Atlas
Model/Module	Stirrer Seal and Guide Maintenance
Author	PP
Date	30/08/2017

Technical Note – Stirrer Seal and Guide Maintenance

The Atlas Stirrer Seal and Guide (Fig 1) can be disassembled without the use of tools for maintenance very easily.



Figure 1

1. Unscrew the lower section of the Stirrer Seal and Guide from the main body (Fig 2)



Figure 2

2. Remove the PEEK section from the Aluminium body (Fig 3)



Figure 3

3. Press down and rotate the central black seal anti-clockwise to release the spring and seal assembly (Fig 4)



Figure 4

4. Remove the central seal and spring assembly (Fig 5)



Figure 5

5. The stirrer seal and guide is now disassembled and ready for seal replacement or general maintenance.